


iINTER-SHOT

<shooting screen>

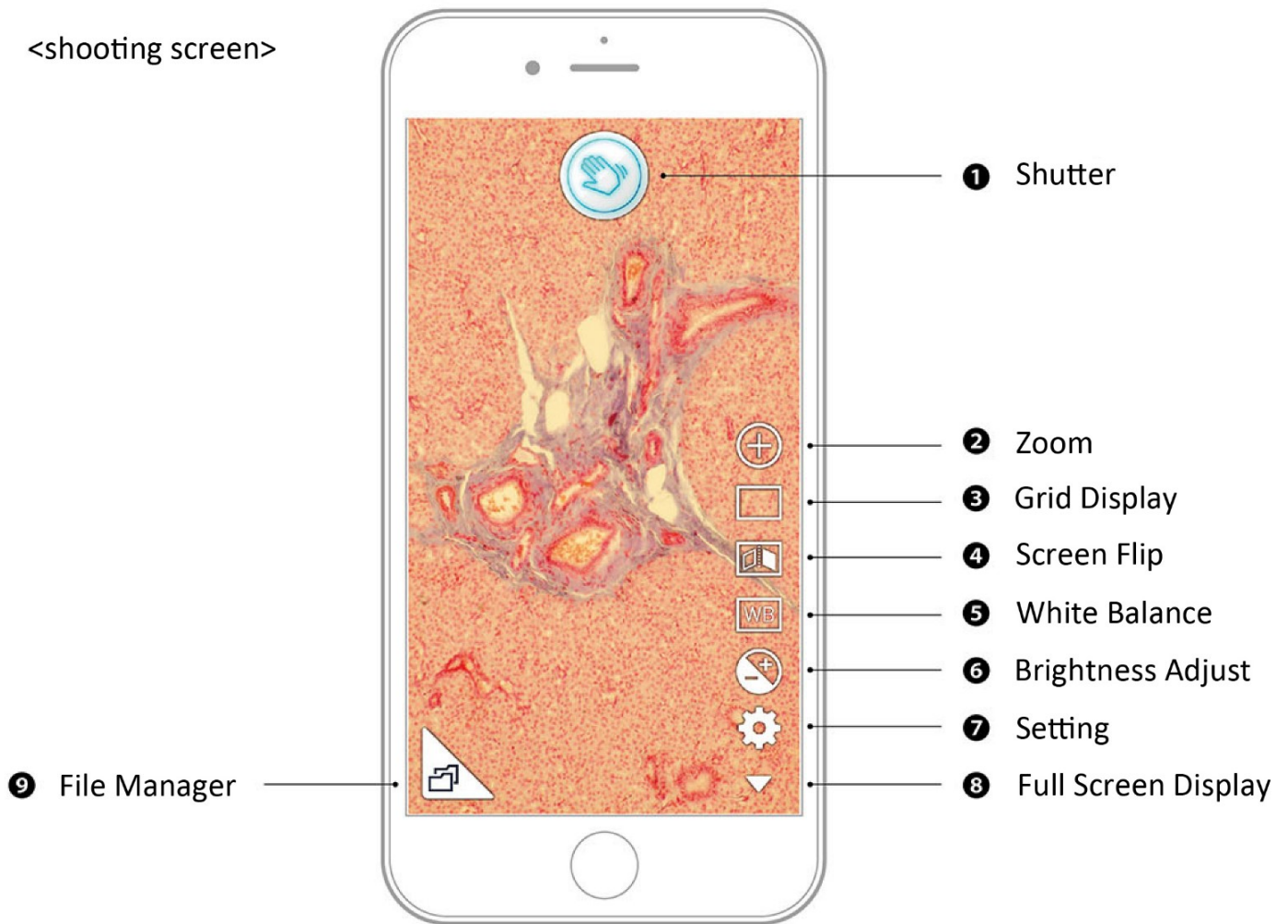

- ① Shutter : Hold hand over the icon for touch-less shutter. Tapping the icon can release the shutter too.
- ② Zoom : Three levels of zooming are available. Convenient for focusing.
- ③ Grid Display : Three types of grid displays are available.
- ④ Screen Flip : Flip screen up side down for microscope which renders reverted image.
- ⑤ White Balance : Manual adjustment with bar gauge. Tap **AUTO** for auto adjustment.
- ⑥ Brightness Adjustment : ISO Sensitivity can be adjusted. Tap **AUTO** for auto adjustment.
- ⑦ Setting : Move to setting menu.
- ⑧ Full Screen Display : Remove all icons for clear full screen view.
- ⑨ File Manager : Move to file manager.

<confirmation screen>

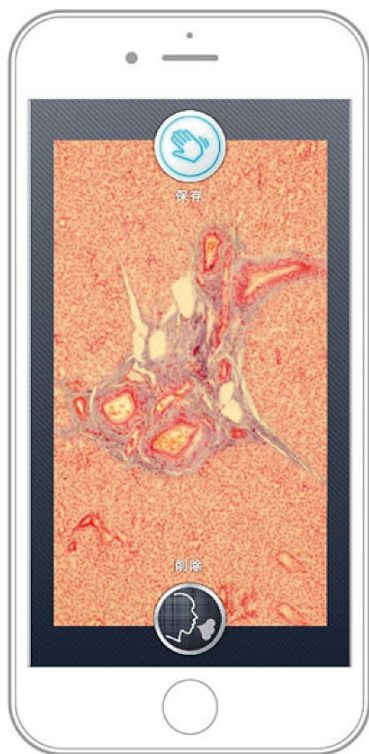


Image will be saved with holdind hand over hand icon after releasing shutter

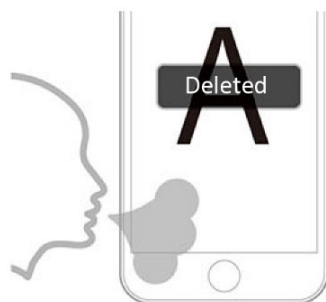


Image will be deleted with breathing on image after releasing shutter

** Breathe delete may disfunction when there is noise in the surrounding. In such case, set it off from setting screen.